# I. JuLI Stage Viewer codec Installation

- Insert the supplied USB drive or go to Dropbox and copy the JuLI Stage Viewer SW' folder on the desktop.
- 2. Double click 'Xvid-1.3.3-20140407' file to install.



3. Select language and press 'OK'.

X Language Selection	
Please select the in	nstallation language
English - English	•
ОК	Cancel

4. Press 'Next' button.



5. Accept the agreement and press 'Next' button.

🔀 Setup		
License Agreement	×	
Please read the following Lic agreement before continuing	ense Agreement. You must accept the terms of this g with the installation.	
GNU	J GENERAL PUBLIC LICENSE	
	Version 2, June 1991	
Copyright (C) 1989, 1991 Free Software Foundation, Inc. 59 Temple Place, Suite 330, Boston, MA 02111-1307 USA Everyone is permitted to copy and distribute verbatim copies of this license document, but changing it is not allowed.		
	Preamble	
Do you accept this license?	<ul> <li>I accept the agreement</li> <li>I do not accept the agreement</li> </ul>	
BitRock Installer	,	
	< Back Next > Cancel	

6. Set installation directory and press 'Next' button.

Setup	
Installation Directory	×
Please specify the directory where Xvid Video Codec will be installed.	
Installation Directory C#Program Files (x86)#Xvid	- <u>6</u>
BitRock Installer	
< Back Next >	Cancel

7. Select update setting and press 'Next' button.



 Select 'DIVX' type essentially and press 'Next' button.

🔀 Setup			
Decoder Configuration			X
Use Xvid to play back also the f	ollowing media types:		
☑ DIVX			
3IVX			
Other MPEG-4			
BitRock Installer			
	< Back	Next >	Cancel



### NanoEnTek Inc. (HQ)

12F, 5, Digital-ro 26-gil, Guro-gu, Seoul, 152-740, Korea Tel : +82-2-6220-7940, Fax : +82-2-6220-7721

#### NanoEnTek USA Inc.

i627 Stoneridge Drive Suite 304, Pleasanton, CA 94588, USA iel : +1-925-225-0108, Fax : +1-925-225-0109

9. **Press 'Next' button** again, then the installation will be started.

🕻 Setup			
Ready to Install			X
Setup is now ready to begin installing Xvid	Video Codec o	n your compu	uter.

10. Xvid Video Codec is being installed.



11. **Press 'Finish' button** to complete codec installation.





### NanoEnTek Inc. (HQ)

12F, 5, Digital-ro 26-gil, Guro-gu, Seoul, 152-740, Korea Tel : +82-2-6220-7940, Fax : +82-2-6220-7721

#### NanoEnTek USA Inc.

5627 Stoneridge Drive Suite 304, Pleasanton, CA 94588, USA Tel : +1-925-225-0108, Fax : +1-925-225-0109

# **II. JuLI Stage Viewer program Installation**

 Make sure the 'JuLI Stage Projects' folder is on the desktop.



- Note: This project data should be considered as a sample data for JuLI Stage Viewer program.
- Double click 'Setup JuLIStage Viewer V.X.X.X.X' file to install.



## 3. Press 'Next' button.



4. Select installation folder and press 'Next' button.

JuLIStage Viewer
Select Installation Folder
The installer will install JuLIStage Viewer to the following folder.
To install in this folder, click "Next". To install to a different folder, enter it below or click "Browse".
<u>F</u> older:  C:₩Program Files (x86)₩NanoEnTek₩JuLIStage Viewer∜   Disk Cost
Install JuLIStage Viewer for yourself, or for anyone who uses this computer:
C Everyone
Just me
Cancel < Back Next >

5. Press 'Next' button again to confirm installation.

Boldmistage viewer	
Confirm Installation	5
The installer is ready to install JuLIStage Viewer on your computer.	
Click "Next" to start the installation.	
Cancel < Back	Next >

# 6. JuLI Stage Viewer is being installed.

🛃 JuLIStage Viewer		
Installing JuLIStage Vie	er	
JuLIStage Viewer is being installed.		
Please wait		
	Cancel < Bac	k Next>
	Cancel < Bac	ik Next >

7. Press 'Close' button to complete the installation.



# Nano**EnTek** sales@nanoentek.com www.nanoentek.com

### NanoEnTek Inc. (HQ)

12F, 5, Digital-ro 26-gil, Guro-gu, Seoul, 152-740, Korea Tel : +82-2-6220-7940, Fax : +82-2-6220-7721

### NanoEnTek USA Inc.

5627 Stoneridge Drive Suite 304, Pleasanton, CA 94588, USA Tel : +1-925-225-0108, Fax : +1-925-225-0109

8. Double click 'JuLIStage Viewer' file on the

desktop to run the JuLI Stage Viewer program..



Select a data directory path. Click 'Browse...'
 button to choose a folder containing projects that you want to import. Choose 'JuLI Stage Projects' folder on the desktop that you saved at the first step. Then press 'OK' button.



 Note: If you have other project folder from the JuLI Stage, copy that project folder in the 'JuLI Stage Projects' folder.
 Note: You have to select folder which contains project folders. Do not select a single project folder.



### NanoEnTek Inc. (HQ)

12F, 5, Digital-ro 26-gil, Guro-gu, Seoul, 152-740, Korea Tel : +82-2-6220-7940, Fax : +82-2-6220-7721

#### NanoEnTek USA Inc.

5627 Stoneridge Drive Suite 304, Pleasanton, CA 94588, USA Tel : +1-925-225-0108, Fax : +1-925-225-0109