Nano EnTek

JuLI Stage Update Instruction

(Offline, Manual update)

1. Open 'JuLI Stage Update' folder.



2. Open 'DirectX Setup' file and double click 'DXSETUP' icon.



3. Select 'I accept the agreement' and click 'Next'.



4. Keep clicking the 'Next' button, then installation will be completed. Click 'Finish'.



5. Double click 'JuLI Stage' icon on the desktop.



6. Go to 'Settings' menu and **click 'Update'** next to the 'Software version'.



7. Select 'JuLI Stage' icon which you want to update and click 'Open' button.

-	
JuLI Stage	<u>O</u> pen



NanoEnTek Inc. (HQ)

12F, 5, Digital-ro 26-gil, Guro-gu, Seoul, 152-740, Korea Tel : +82-2-6220-7940, Fax : +82-2-6220-7721

NanoEnTek USA Inc.

5627 Stoneridge Drive Suite 304, Pleasanton, CA 94588, USA Tel : +1-925-225-0108, Fax : +1-925-225-0109

Nano**EnTek**

8. Click 'OK'. Then it will reboot the JuLI Stage.



9. Click 'OK' and it will be rebooted to install.



10. Click 'OK' again.



11. Go to 'Settings' menu and **click 'Update'** next to the 'Firmware version'.



anoEnle

sales@nanoentek.com www.nanoentek.com **12.** It updates and reboots automatically.



13. The **updated version** will be displayed.

Version Auto ch	eck V	
Software information		Prev
Software Version	1.0.2.0 Update	/iew
New Version	Unknown Check Update	
Firmware information		
Firmware Version	H 1.00 F 0.13 Update	•
New Version	Unknown Check Update	ption
Objective Lens	⊙ x4	
Data Path	E:\JuLIStage Set	
Total Space Free Space Used Space	471,250,940 MByte 470,636,540 MByte 00.13 %	Data
Mouse control	Type 1 L mouse : Select R mouse : Move L mouse : Select	Sett
Serial	JS141203-008	tings
Admin	Admin Option Locking Set	_ ×

NanoEnTek Inc. (HQ)

12F, 5, Digital-ro 26-gil, Guro-gu, Seoul, 152-740, Korea Tel : +82-2-6220-7940, Fax : +82-2-6220-7721

NanoEnTek USA Inc.

5627 Stoneridge Drive Suite 304, Pleasanton, CA 94588, USA Tel : +1-925-225-0108, Fax : +1-925-225-0109